[www.ricerca.it](http://www.ricerca.it)

**PROMOZIONE**

Ver. 2024/02

**VALIDITA'**

dal 01/08/2024 al 31/12/2024

**CODICE**

FBSMinis\_2024

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### The Essentials at a Glance

	FBS STANDARD	FBS MINIS
<b>Aliquoting</b>	Yes	No
<b>Working time</b>	Thawing bottle: 30 min Aliquoting: 15 min Water bath: 15 min	Water bath: 15 min
<b>Plastic waste</b>	100 %	70 %
<b>Freeze-thaw cycles</b>	2-3	1
<b>Labeling</b>	Manually & error-prone	Distinct label with Cat. No., Lot & Expiry date

Codice	Descrizione	Conf.	Listino	Promozione
<b>R110-FBS-12F</b>	Fetal Bovine Serum, Collected in South America	10x50 ml	265,00€	212,00€
<b>R110-FBS-HI-12F</b>	Fetal Bovine Serum (FBS Minis), Heat Inactivated, Collected in South America	10x50 ml	276,00€	220,80€
<b>R110-FBS-11F</b>	Fetal Bovine Serum Advanced, Collected in South America	10x50 ml	233,00€	186,40€
<b>R110-FBS-HI-11F</b>	Fetal Bovine Serum Advanced (FBS Minis), Heat Inactivated, Collected in South America	10x50 ml	239,00€	191,20€
<b>R110-FBS-16F</b>	FBS Xtra, Fetal Bovine Serum, Collected in South America	10x50 ml	217,00€	173,60€
<b>R110-FBS-HI-16F</b>	FBS Xtra, Fetal Bovine Serum (FBS Minis), Heat Inactivated, Collected in South America	10x50 ml	220,00€	176,00€

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## Frequently Asked Questions

### What is the shelf life of FBS Minis?

Our FBS Minis have a shelf life of 5 years from the manufacturing date of the original batch.

However, please note that there may be differences in shelf life depending on the specific variant.

### Can I heat-inactivate the FBS Minis?

Yes, the FBS Minis can be heat-inactivated without any issues, and as usual.

Additionally, we offer the option to provide you with heat-inactivated FBS Minis, offering a further time-saving advantage.

### How can the FBS Minis be thawed?

Your serum can be thawed overnight at +4 °C to +8 °C (controlled conditions) or in a water bath at +37 °C while gently swirling occasionally.

Once the serum is fully thawed, gently invert the bottle to ensure the mixing of all components. Carefully add the serum to your culture medium or the flask to be used.

### How do I secure my bottles in the water bath?

Adjust your water level to accommodate the bottles without extra attachments. Thus, you do not need any other supporting attachments for your bottles. If this is not sufficient, you can weigh down your Minis with common utensils. Make sure that the method you choose provides a stable hold for the bottles.

Moreover, we recommend a controlled thawing process at +4 °C to +8 °C.

### What material are the bottles made from?

Our FBS Minis bottles are constructed from modified polyethylene terephthalate with glycol (PETG). They are specifically designed for medical use and allow easy sterilization. The material is characterized by its recyclability, which makes PETG an extremely eco-friendly option.

### Can I transfer my Minis by pouring, or do I HAVE TO use a pipette?

Yes, pouring your Minis is a viable option, which at the same time contributes to the reduction of plastic waste. However, it is advisable to be careful during the pouring process and transfer the contents precisely to avoid spilling.

In particular, our Minis are well-suited for preparing medium with 10% FBS since you typically use the entire bottle, preventing any backflow or contamination risks.

### How should I label my aliquots?

Do not worry about using pens or inconvenient stickers that might fall off or smudge. Each of our Minis already comes with a label showing you all the important information.

In addition, the storage box provides the flexibility to add other relevant data, such as the user or workgroup name, for your convenience.

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